
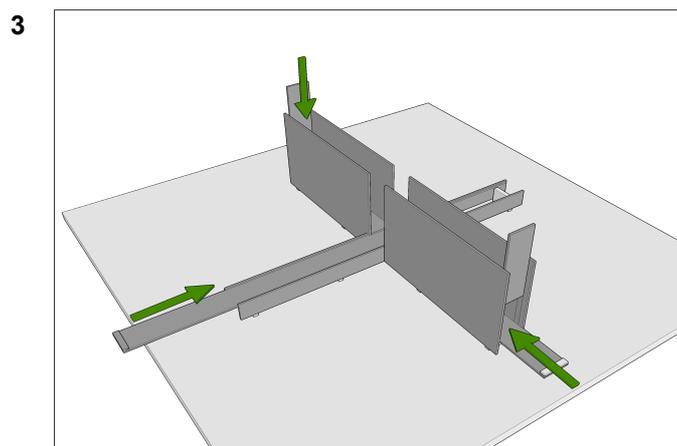
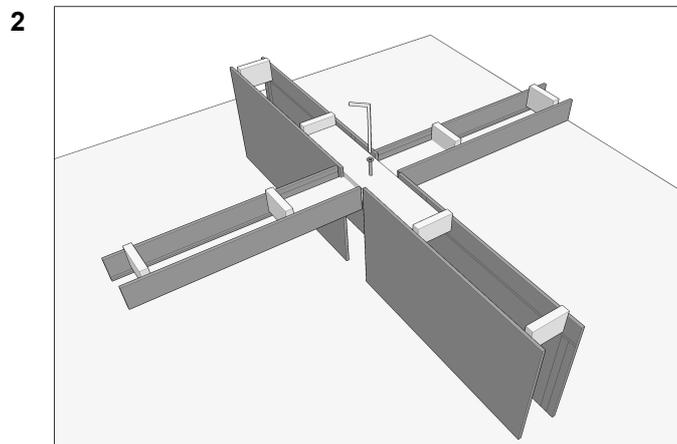


Assembly Instructions

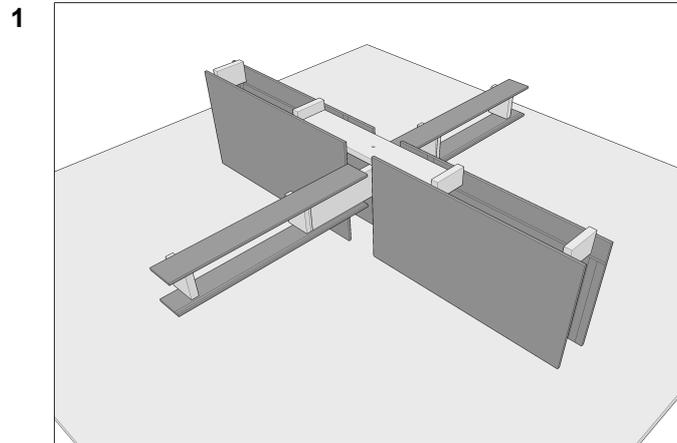

Assembling instructions

Take away the protective film from all details.

If the cross-piece of the cart has been shipped non-assembled, please assemble it using four screws and four nuts as it is shown in the picture below.

Attach a leg with a wheel to every end of a cross-piece. Every leg is attached by means of three screws and three nuts with a nylon ring.

Connect the opposite legs with long straight pipes using the type ❶ fasteners. Do not tighten the screws of the type ❶ fasteners, allow them to move freely along the legs.

Attach the type ❷ fastener to the crossing of straight pipes. Do not tighten the screw.

Put the cart on a flat surface and ensure that all wheels are touching it even when rotating the cart.

Tighten the type ❷ fastener.

Position both pipes and their type ❶ fasteners 3-5cm above the wheels (the lower pair of them). Tighten the type ❶ fasteners.

If the cart was assembled correctly but it is still unsteady on an uneven surface just rotate the cart until it becomes steady.

